

Double-Strand Specific DNase (dsDNase)

Product Data

Alternative names

Shrimp DNase

Source

Produced in a *Pichia pastoris* strain expressing a recombinant gene for a nuclease from *Pandalus borealis* (arctic shrimp).

Concentration

1-50 U/ μ l

Specific activity

Ca. 475 000 Unit/mg

Unit definition

One unit increases the absorbance at 260 nm by 0.001 OD per min at 25°C and pH 5.0 with large molecular weight DNA as the substrate according to the assay method of Kunitz.

Storage

Store at -20°C

Stable at 4°C, and can also be stored for prolonged time in a refrigerator.

Storage buffer

20 mM Tris-HCl (pH 7.5 at 25°C)

2.0 mM MgCl₂

10 mM NaCl

0.01% (v/v) Triton X-100

50% (v/v) glycerol

Protocol recommendations

For preferential degradation of dsDNA and no RNase activity, use a reaction buffer containing 20 mM Tris-HCl (pH 8.0) and 2.5 mM MgCl₂.

Activators

Mg²⁺ ions (up to 10 mM) are required for maximum activity.

Optimum pH: 7.5

Optimum reaction temperature: 40°C

Purity

The enzyme is purified to apparent homogeneity by SDS-PAGE.

Properties

dsDNase is an endonuclease that cleaves phosphodiester linkages in DNA to yield oligonucleotides with 5'-phosphate and 3'-hydroxyl termini. dsDNase has a very high specific activity, estimated 30 times higher than bovine DNase I, and it is heat labile. dsDNase has a particularly strong preference for double-stranded DNA (dsDNA). In the presence of magnesium as only divalent cation and using oligos as a substrate; the activity towards dsDNA is 5000-fold higher than towards ssDNA. The enzyme can therefore be used to specifically degrade dsDNA, leaving ssDNA essentially intact.

Applications

Removal of contaminating DNA before PCR or RT-PCR leaving single-stranded primers and RNA intact.

Heat inactivation

The enzyme is easily heat-inactivated at 65°C for 15 min in the presence of 1 mM DTT.

For further information contact:

Gerd Nilsen

gn@arcticzymes.com

+47 45 47 00 77

Disclaimer Certain applications of ArcticZymes AS products may require licenses from others. It is the expressed duty of any receiver of ArcticZymes AS products to acquire such licenses, if necessary. In no event shall ArcticZymes AS be liable for claims for any damages, whether direct, incidental, foreseeable, consequential, or special (including but not limited to loss of use, revenue or profit), arising due to the violation of third parties Intellectual Property Rights by any receiver of ArcticZymes AS products. ArcticZymes AS products may be covered by pending or issued patents, designs or design applications and/or trademarks or trademark applications or any other registered or unregistered Intellectual Property Right.

